

Iron Low Range & Manganese Low Range Photometer





Dear
Customer,Thank you for choosing a Hanna Instruments product.Please read this instruction manual carefully before using this instrument.
This manual will provide you with the necessary information for correct
use of this instrument, as well as a precise idea of its versatility.
If you need additional technical information, do not hesitate to e-mail
us at tech@hannainst.com or view our worldwide contact list at
www.hannainst.com.

All rights are reserved. Reproduction in whole or in part is prohibited without the written consent of the copyright owner, Hanna Instruments Inc., Woonsocket, Rhode Island, 02895, USA.

TABLE OF CONTENTS

1. PRELIMINARY EXAMINATION	
2. SAFETY MEASURES	
3. ABBREVIATIONS	5
4. SPECIFICATIONS	
5. DESCRIPTION	7
5.1. GENERAL DESCRIPTION & INTENDED USE	7
5.2. FUNCTIONAL DESCRIPTION	8
5.3. PRECISION & ACCURACY	9
5.4. PRINCIPLE OF OPERATION	9
5.5. OPTICAL SYSTEM	10
6. GENERAL OPERATIONS	
6.1. METER VALIDATION: CAL CHECK & CALIBRATION	11
6.2. CHEMICAL FORMULA & UNIT CONVERSION	13
6.3. GLP	
6.4. LOGGING DATA & LOG RECALL	13
6.5. GENERAL SETUP	
6.6. REAGENTS & ACCESSORIES	
6.7. CONTEXTUAL HELP	17
6.8. BATTERY MANAGEMENT	18
7. PHOTOMETER	
7.1. METHOD SELECTION	
7.2. COLLECTING & MEASURING SAMPLES AND REAGENTS	
7.3. CYLINDER & CUVETTE PREPARATION	
8. METHOD PROCEDURE	
8.1. IRON LOW RANGE	
8.2. MANGANESE LOW RANGE	
9. WARNING & ERROR DESCRIPTIONS	
10. BATTERY REPLACEMENT	
11. ACCESSORIES	
11.1. REAGENT SETS	
11.2. OTHER ACCESSORIES	
CERTIFICATION	
RECOMMENDATIONS FOR USERS	
WARRANTY	34

1. PRELIMINARY EXAMINATION

Remove the instrument and accessories from the packaging and examine it carefully. For further assistance, please contact your local Hanna Instruments Office or email us at tech@hannainst.com.

Each H197742C is delivered in a rugged carrying case and is supplied with:

- Sample cuvette (2 pcs.)
- Sample cuvette cap (2 pcs.)
- Plastic stopper (2 pcs.)
- A ZERO CAL Check Cuvette A
- H197742B CAL Check Cuvette B for Iron Low Range
- HI97748B CAL Check Cuvette B for Manganese Low Range
- Cloth for wiping cuvettes
- Scissors
- 1.5V AA Alkaline batteries
- CAL Check standard certificate
- Instrument quality certificate
- Instruction manual

Each H197742 is delivered in a cardboard box and is supplied with:

- Sample cuvette (2 pcs.)
- Sample cuvette cap (2 pcs.)
- Plastic stopper (2 pcs.)
- 1.5V AA Alkaline batteries
- Instrument quality certificate
- Instruction manual

Note: Save all packing material until you are sure that the instrument works correctly. Any damaged or defective item must be returned in its original packing material with the supplied accessories.

2. SAFETY MEASURES



- The chemicals contained in the reagent kits may be hazardous if improperly handled.
- Read the Safety Data Sheets (SDS) before performing tests.
- Safety equipment: Wear suitable eye protection and clothing when required, and follow instructions carefully.
- Reagent spills: If a reagent spill occurs, wipe up immediately and rinse with plenty of water. If reagent contacts skin, rinse the affected area thoroughly with water. Avoid breathing released vapors.
- Waste disposal: For proper disposal of reagent kits and reacted samples, contact a licensed waste disposal provider.

3. ABBREVIATIONS

μ g/L	micrograms per liter (ppb)
mg/L	milligrams per liter (ppm)
mL	milliliter
°C	degree Celsius
°F	degree Fahrenheit
EPA	US Environmental Protection Agency
GLP	Good Laboratory Practice
HDPE	High Density Polyethylene
LED	Light Emitting Diode
LR	Low Range
NIST	National Institute of Standards and Technology

4. SPECIFICATIONS

	Range	0.00 to 1.60 mg/L (as Fe)	
Iron LR	Resolution	0.01 mg/L	
Ifon LK	Accuracy	\pm 0.01 mg/L \pm 8% of reading at 25 °C	
	Method	Adaptation of the TPTZ Method	
	Range	0 to 300 µg/L (as Mn)	
	Resolution	1 μg/L	
Manganese LR	Accuracy	\pm 10 μ g/L \pm 3% of reading at 25 °C	
	Method	Adaptation of the 1-(2-Pyridylazo)-2-Naphthol PAN Method	
	Light source	Light Emitting Diode	
	Bandpass filter	575 nm	
Measurement	Bandpass filter bandwidth	8 nm	
System	Bandpass filter wavelength accuracy	±1.0 nm	
	Light detector	Silicon photocell	
Cuvette type		Round 24.6 mm diameter (22 mm inside)	
	Auto logging	50 readings	
	Display	128 x 64 pixel B/W LCD with backlight	
	Auto-off	After 15 minutes of inactivity (30 minutes before a READ measurement)	
	Battery type	1.5 V AA Alkaline (3 pcs.)	
Additional	Battery life	> 800 measurements (without backlight)	
Specifications	Environment	0 to 50 °C (32 to 122 °F); 0 to 100% RH, non-serviceable	
	Dimensions	142.5 x 102.5 x 50.5 mm (5.6 x 4.0 x 2.0")	
	Weight (with batteries)	380 g (13.4 oz.)	

5. DESCRIPTION

5.1. GENERAL DESCRIPTION & INTENDED USE

The HI97742 is an auto-diagnostic portable photometer that benefits from Hanna's years of experience as a manufacturer of analytical instruments. It has an advanced optical system that uses a Light Emitting Diode (LED) and a narrow band interference filter that allows for accurate and repeatable readings.

The optical system is sealed from outside dust, dirt and water. The meter uses an exclusive positive-locking system to ensure that the cuvettes are placed into the holder in the same position every time.

With the CAL Check[™] functionality, users are able to validate the performance of the instrument at any time and apply a user calibration (if necessary). Hanna Instruments CAL Check cuvettes are made with NIST traceable standards.

The built-in tutorial mode guides users step-by-step through the measurement process. It includes all steps required for sample preparation, the required reagents and quantities.


The HI97742 meter measures the iron content in water samples from 0.00 to 1.60 mg/L (ppm) and manganese content in water samples in the 0 to 300 μ g/L (ppb).


The method for iron is an adaptation of the TPTZ Method. The method for manganese is an adaptation of the 1-(2-Pyridylazo)-2-Naphthol PAN Method.

The H197742 photometer is a compact and versatile meter suitable for field or bench measurements, featuring a:

- Sophisticated optical system
- Meter validation using certified CAL Check cuvettes
- Tutorial mode guides the user step-by-step
- Auto logging
- Waterproof IP67, floating case
- GLP features

5.2. FUNCTIONAL DESCRIPTION

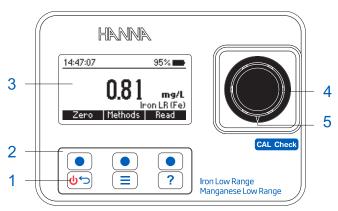


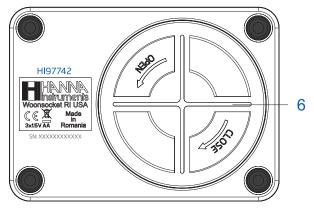

1. ON/OFF power button3. Liquid Crystal Display (LCD)5. Indexing mark2. Keypad4. Cuvette holder6. Battery cover

Keypad Description

The keypad contains 3 direct keys and 3 functional keys with the following functions:

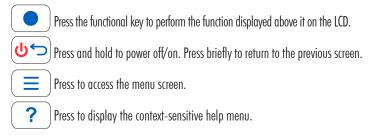


5.3. PRECISION & ACCURACY

Precision is how closely repeated measurements are to one another. Precision is usually expressed as standard deviation (SD).

Accuracy is defined as the closeness of a test result to the true value and is method specific.

Although good precision suggests good accuracy, precise results can be inaccurate. The figure explains these definitions.

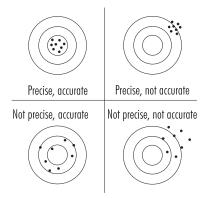

5.4. PRINCIPLE OF OPERATION


Absorption of light is a typical phenomenon of interaction between electromagnetic radiation and matter. When a light beam crosses a substance, some of the radiation may be absorbed by atoms, molecules or crystal lattices. Photometric chemical analysis is based on specific chemical reactions between a sample and reagent to produce a light-absorbing compound.

If pure absorption occurs, the fraction of light absorbed depends both on the optical path length through the matter and on the physical-chemical characteristics of the substance according to the Lambert-Beer Law. If all other factors are constant, the concentration "c" can be calculated form the absorbance of the substance.

$$\begin{array}{c} \text{-log I/I}_{\mathrm{o}} = \epsilon_{\lambda} \, \text{c} \, \text{d} \\ & \text{or} \\ \text{A} = \epsilon_{\lambda} \, \text{c} \, \text{d} \end{array}$$

- I_{o} = intensity of incident light beam
- I = intensity of light beam after absorption
- ϵ_{λ} = molar extinction coefficient at wavelength λ
- c = molar concentration of the substance
- d = optical path through the substance

5.5. OPTICAL SYSTEM

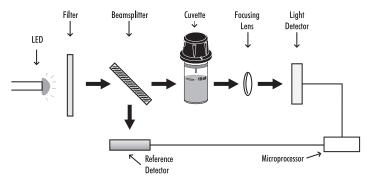

Instrument Block Diagram

The internal reference system (reference detector) of the H197742 photometer compensates for any drifts due to power fluctuations or ambient temperature changes, providing a stable source of light for your blank (zero) measurement and sample measurement.

LED light sources offer superior performance compared to tungsten lamps. LEDs have a much higher luminous efficiency, providing more light while using less power. They also produce little heat, which could otherwise affect electronic stability. LEDs are available in a wide array of wavelengths, whereas tungsten lamps have poor blue/ violet light output.

Improved optical filters ensure greater wavelength accuracy and allow a brighter, stronger signal to be received. The end result is higher measurement stability and less wavelength error.

A focusing lens collects all of the light that exits the cuvette, eliminating errors from cuvette imperfections and scratches, eliminating the need to index the cuvette.

6. GENERAL OPERATIONS

6.1. METER VALIDATION: CAL CHECK & CALIBRATION

Validation of the H197742 involves verifying the concentration of the certified CAL Check standards. The CAL Check screen guides the user step-by-step through the validation process and user calibration (if necessary).

WARNING: Do not use any solutions or standards other than the Hanna Instruments[®] CAL Check Standards. For accurate validation and calibration results, please perform these at room temperature, 18 to 25 °C (64.5 to 77.0 °F).

Note: CAL Check Standards will not read the specified value in measurement mode. Protect the CAL Check cuvettes from direct sunlight by keeping them in the original packing. Store between 5 and 30 °C (41 to 86 °F), do not freeze.

To perform a CAL Check:



The "Not Available" message or the date, time and status of the last CAL Check will be displayed on the screen.



Note: CAL Check & Calibration is for the selected method.

- 2. Press **Check** to start a new CAL Check. Press the 🕑 ⊃ key at any time to abort the validation process.
- 3. Use the functional keys to enter the certificate value of the calibration standard found on the CAL Check Standard Certificate. Press **Next** to continue.



Note: This value will be saved in the instrument for future validation. If a new set of calibration standards is obtained please update the certificate value.

5.

 Insert the A ZERO CAL Check Cuvette A then press Next to continue. The "Please wait..." message will be displayed during the measurement.

95% 💼




Insert the CAL Check Cuvette B for the selected method (HI97742B for Iron LR or
H197748B for Manganese LR) then press Next to continue. The "Please wait"
message will be displayed during the measurement.



- 6. When the CAL Check is complete the display will show one of the following messages and the value obtained during the measurement:
 - "PASSED": The measured value is within the accuracy specification, no user calibration is required.



 "OUT OF SPECIFICATION" and Calibrate is available: The measured value is near the expected value. To update the user calibration press Calibrate. Press Accept to confirm or Cancel to return to the previous screen.

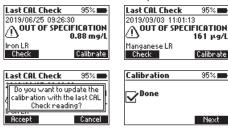

 "OUT OF SPECIFICATION": A user calibration is not allowed, the measured value is outside of the tolerance window. Check the certified value, expiration date and clean the outside of the cuvette. Repeat the CAL Check procedure. If this error continues contact your nearest Hanna Instruments Customer Service Center.



6.2. CHEMICAL FORMULA & UNIT CONVERSION

Chemical formula and unit conversion factors are pre-programmed into the instrument and are method specific. In order to view the displayed result in the desired chemical formula, for the manganese LR method, enter menu by pressing and use the functional keys to select *Chemical Form*. Press **Select** to change the displayed chemical formula. Use the functional keys to highlight the desired chemical formula and press **Select**. The selected formula will be saved when the instrument is powered off.



6.3. GLP

Press the \equiv key to enter the menu. Use the functional keys to select *GLP* and press **Select**. Good Laboratory Practice (GLP) shows the date and time of the last user calibration (if available) or factory calibration. To erase the last user calibration and to clear the CAL Check press **Clear** and follow the prompts. Press **Yes** to erase and return to the factory calibration data or **No** to exit the clear procedure.


Last Calibration	95% 💼
2019/10/11 10:23:21	
Iron LR	
	Clear

6.4. LOGGING DATA & LOG RECALL

The instrument features a data autolog function to help users keep track of all measurements. Every time a measurement is made the data is automatically saved. The data log can hold 50 individual measurements. When the data log is full (50 data points) the meter will rewrite the oldest data point.

Viewing and deleting the data is possible using the Log Recall menu.

Press the (\equiv) key to enter the menu. Use the functional keys to select *Log Recall* and press **Select**.

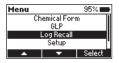

Use the functional keys to highlight a log and press **Info** to view additional information about the log. From this screen **Next** and **Previous** can be used to view other logs.

og Reca	all 4/12	95% 💼	Log Recall 95%
10/17	0.33 mg/L	Fe .	0.12 mg/L Fe
10/17	0.12 mg/L	Fe	Iron LR
10/17	220 µg/L	Mn	2019/10/17 13:05:22
10/17	57 µg/L	Mn	
	I ▼ 1	Info	Delete Previous Next

Press **Delete** to erase logged data. After pressing **Delete** a prompt on display is asking for confirmation.



Press No or the 0 key to return to the previous screen.

Press Yes to delete selected log.

Press **Del All** to erase all the logged data. If **Del All** is pressed follow the prompt to confirm. Press **Yes** to delete all logged data, **No** or the **O** so the **O** so the **D** several.

6.5. GENERAL SETUP

Press the 📃 key to enter the menu. Use the functional keys to select *Setup* and press **Select**. Use the functional keys to highlight desired option.

Backlight

Option: 0 to 100 %


Press **Modify** to access the backlight intensity. Use the functional keys to increase or decrease the value. Press **Accept** to confirm or the key to return to the *Setup* menu without saving the new value.

Setup	95% 💼	Backlight	95%
Backlight	50%	0%	10
Contrast	50%	0%	10
Date / Time	07:23:32	503	%
Time Format	24-hour		
▲ ▼	Modify		Accep

Contrast

Option: 0 to 100 %

Press **Modify** to change the display's contrast. Use the functional keys to increase or decrease the value. Press **Accept** to confirm the value or the *Setup* menu without saving the new value.

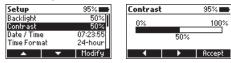

Date & Time

Press **Modify** to change the date and time. Press the functional keys to highlight the value to be modified (year, month, day, hour, minute or second). Press **Edit** to modify the highlighted value. Use the functional keys to change the value.

Press Accept to confirm or the 0 key to return to the previous screen.



Time Format

Option: AM/PM or 24-hour

Press the functional key to select the desired time format.

Setup	95% 💼
Contrast	50%
Date / Time	07:25:14
Time Format	24-hour
Date Format	YYYY/MM/DD
▲	👻 🛛 AM/PM

Date Format

Press **Modify** to change the date format. Use the functional keys to select the desired format. Press **Accept** to confirm or the key to return to the *Setup* menu without saving the new format.

95%1

at

Setup		95% 💼	Date For	
Date / Time	2	07:25:26	DD/MM/YY	
Time Form		24-hour	MM/DD/YY	
Date Form		Y/MM/DD	YYYY/MM	
Decimal Se	parator	•	YYYY-MM	-D
	•	Modify	•	

Decimal Separator

Option: Comma (,) or Period (.)

Press the functional key to select the desired decimal separator. The decimal separator is used on the measurement screen.

Setup		95% 💼
Time Form	at 24-hour	
Date Form		
Decimal Se	parator 🔹	
Language		English
▲	•	,

Language

Press **Modify** to change the language. Use the functional keys to select the desired language. Press **Accept** to choose one of the languages installed.



Beeper

Option: Enable or Disable

When enabled, a short beep is heard every time a key is pressed. A long beep alert sounds when the pressed key is not active or an error is detected. Press the functional key to enable or disable the beeper.

Setup		95% 💼
Decimal Sep	arator	•
Language		English
Beep On		
Tutorial		
	•	Disable

Tutorial

Option: Enable or Disable

When enabled, the user will be guided step-by-step through the measurement procedure.

Setup		95% 💼
Language		English
Beep On		
Tutorial		
Meter Info	rmation	1
	•	Enable

Meter Information

Press **Select** to view the model, serial number, firmware version and selected language. Press the version between the setup menu.

Setup	95% 💼	Meter Infor	mation
Beep On		Model	HI97742
Tutorial		Serial #	901290005111
1eter Information		Firmware	v1.00
Restore factory setti	ngs	Language	English v1.0
▲ ▼	Select	www.hannainst.com	

Restore Factory Settings

Press **Select** to reset to factory settings. Press **Accept** to confirm or **Cancel** to exit without restoring the factory settings.

Setup 95% 🗰	Setup 95%
Beep On 🛛	Beep On 🗖
Tutorial 🛛 🗖	Do you want to restore
Meter Information	default settings?
Restore factory settings	hestore ractory settings
🔺 🔻 Select	Accept Cancel


6.6. REAGENTS & ACCESSORIES

Press the (\equiv) key to enter the menu. Use the functional keys to select *Reagents / Accessories* and press **Select** to access a list of reagents and accessories. To exit press the 0 key.

95% 💼	Accessories	95%
П	** REAGENT SETS **	
all 📗	Iron LR	
	HI93746-01	
essories	Reagents for 50 tests	
Select	-	
	call p cessories	sall w** REAGENT SETS ** Iron LR HI33746-01 Reagents for 50 tests

6.7. CONTEXTUAL HELP

The H197742 offers an interactive contextual help mode that assists the user at any time. To access the help screen press the ? key.

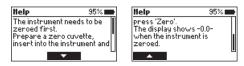


The instrument will display additional information related to the current screen. To read all the available information, scroll the text using the functional keys.

To exit help mode press the 0 or the 2 key and the meter will return to the previous screen.


6.8. BATTERY MANAGEMENT

The meter will perform an auto-diagnostic test when it is powered on. During this test, the Hanna Instruments[®] logo will appear on the LCD. If the auto-diagnostic test was successful, the meter is ready for use. The battery icon on the LCD will indicate the battery status:


• battery full

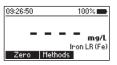

• battery below 10%, replace the batteries soon

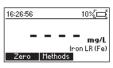

• battery is low, replace the batteries with new ones

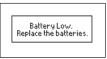

To conserve battery, the meter will turn off automatically after 15 minutes of inactivity. If a zero reading has been done but not a read, auto-off time is increased to 30 minutes.

7. PHOTOMETER

7.1. METHOD SELECTION

Press **Methods** when in measurement mode to access the list of methods. Use the functional keys to highlight the desired method then press **Select**.

Methods	95% 💼
Iron LR Manganese LR	
in any an are an	
▲ I ▼	Select

The selected method will be saved when the instrument is powered off.

7.2. COLLECTING & MEASURING SAMPLES AND REAGENTS


Proper Use of Powder Packet

- 1. Use scissors to open the powder packet.
- 2. Push the edges of the packet to form a spout.
- 3. Pour out the content of the packet.




Proper Use of Syringe

- 1. Push the plunger completely into the syringe and insert the tip into the solution.
- Pull the plunger up until the lower edge of the seal is exactly on the mark for the desired volume.
- Take out the syringe and clean the outside of the syringe tip, be sure that no drops are hanging on the tip of the syringe. Then, keeping the syringe in vertical position above the cuvette, push the plunger down into the syringe, the desired volume has been delivered into the cuvette.

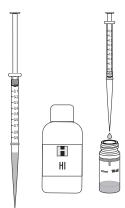

Proper Use of Dropper Bottle

- 1. Tap the dropper on the table several times and wipe the outside of the tip with a cloth.
- 2. Always keep the dropper bottle in a vertical position while dosing the reagent.

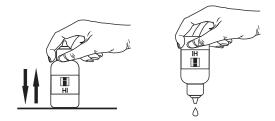

7.3. CYLINDER & CUVETTE PREPARATION

Proper mixing is very important for reproducibility of the measurements. The proper mixing technique is listed in the method procedure.


(a) Invert the cuvette a couple of times or for a specified time: hold the cuvette in the vertical position. Turn the cuvette upside-down and wait for all of the solution to flow to the cap end, then return the cuvette to the upright vertical position and wait for all of the solution to flow to the cuvette bottom. This is one inversion. The correct speed for this mixing technique is 10-15 complete inversions in 30 seconds. This mixing technique is indicated with "invert to mix" and the following icons:



(b) The mixing method is indicated with "shake gently" using one of the following icons:



(c) The mixing method is indicated with "shake vigorously" using one of the following icons:

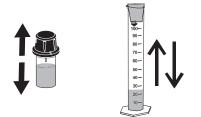

In order to avoid reagent leaking and to obtain more accurate measurements, close the cuvette first with the supplied HDPE plastic stopper _____ and then the black cap.



Whenever the cuvette is placed into the measurement holder, it must be dry outside and free of fingerprints, oil or dirt. Wipe it thoroughly with HI731318 microfiber cleaning cloth or a lint-free wipe prior to insertion.



Shaking the cuvette can generate bubbles in the sample, causing higher readings. To obtain accurate measurements, remove such bubbles by swirling or by gently tapping the cuvette.

Do not let the reacted sample stand too long after reagent has been added. For best accuracy, respect the timings described in the method.

It is possible to take multiple readings in a row, but it is recommended to take a new zero reading for each sample and to use the same cuvette for zeroing and measurement when possible.

Discard the sample immediately after the reading has been taken, or the glass might become permanently stained.

All the reaction times reported in this manual are at 25 °C (77 °F). In general, the reaction time should be increased for temperatures lower than 20 °C (68 °F), and decreased for temperatures higher than 25 °C (77 °F).

8. METHOD PROCEDURE

8.1. IRON LOW RANGE

REQUIRED REAGENTS

Code	Description	Quantity
HI93746-0	Iron LR Reagent	2 packets

REAGENT SETS

HI93746-01 Iron LR Reagent - 50 tests HI93746-03 Iron LR Reagent - 150 tests For other accessories see ACCESSORIES section.


MEASUREMENT PROCEDURE

• Select the Iron LR method using the procedure described in the METHOD SELECTION section.

Note: If tutorial mode is disabled, follow the measurement procedure below. If the tutorial mode is enabled, press **Measure** and follow the messages on the screen.

• Fill one graduated mixing cylinder with deionized water up to the 25 mL mark.


- Add one packet of H193746-0 Iron LR Reagent, close the cylinder with a rubber stopper and shake vigorously for 30 seconds. This is the blank.
- Fill a cuvette with 10 mL of the blank (up to the mark). Replace the plastic stopper and the cap.

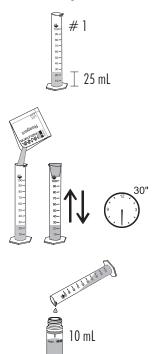

• Insert the cuvette into the holder and ensure that the notch on the cap is positioned securely in the groove.



 Press Zero. The display will show "-0.0-" when the meter is zeroed and ready for measurement.

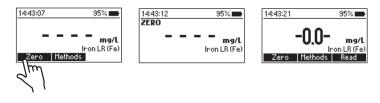


- Remove the cuvette.
- Fill another graduated mixing cylinder with the sample up to the 25 mL mark.



 Add one packet of H193746-0 Iron LR Reagent, close the cylinder with a rubber stopper and shake vigorously for 30 seconds. This is the reacted sample.


• Fill a cuvette with 10 mL of the reacted sample (up to the mark). Replace the plastic stopper and the cap.

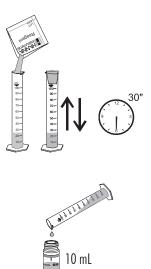


 Insert the cuvette into the holder and ensure that the notch on the cap is positioned securely in the groove.



 Press Read and the display will show a 30 second countdown prior to the measurement. To skip the timer press Read twice. When the timer ends the meter will perform the reading. The instrument displays concentration in mg/L of iron (Fe).

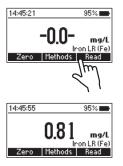



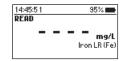

INTERFERENCES

Interference may be caused by:

- Manganese above 50.0 mg/L
- Cadmium, Molybdenum above 4.0 mg/L
- Cyanide above 2.8 mg/L
- Chromium(VI) above 1.2 mg/L
- Nickel above 1.0 mg/L
- Nitrite ion above 0.8 mg/L
- Copper above 0.6 mg/L
- Mercury above 0.4 mg/L
- Chromium(III) above 0.25 mg/L
- Cobalt above 0.05 mg/L
- Sample pH should be between 3 and 4 to avoid fading or turbidity formation

8.2. MANGANESE LOW RANGE

REQUIRED REAGENTS

Code	Description	Quantity
HI93748A-0	Manganese LR Reagent A	2 packets
HI93748B-0	Manganese LR Reagent B	0.40 mL
HI93748C-0	Manganese LR Reagent C	2 mL
HI93703-51	Dispersing Agent	6 drops

REAGENT SETS


HI93748-01 Manganese LR Reagent - 50 tests HI93748-03 Manganese LR Reagent - 150 tests For other accessories see ACCESSORIES section.

MEASUREMENT PROCEDURE

• Select the Manganese LR method using the procedure described in the METHOD SELECTION section.

Note: If tutorial mode is disabled, follow the measurement procedure below. If the tutorial mode is enabled, press **Measure** and follow the messages on the screen.

• Fill one cuvette (#1) with 10 mL of deionized water (up to the mark).

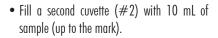

 Add one packet of H193748A-O Manganese LR Reagent A to each cuvette. Replace the plastic stoppers and the caps. Shake gently until completely dissolved.







 Add 0.2 mL of the H193748B-0 Manganese LR Reagent B to each cuvette. Replace the plastic stoppers and the caps. Invert gently to mix for about 30 seconds.

- Add 1 mL of the H193748C-0 Manganese LR Reagent C to each cuvette, replace the plastic stoppers and the caps. Shake gently for 15 seconds.

1 mL

0.2 mL


HI93748B-0

Agent to each cuvette. Replace the plastic stoppers and the caps. Invert gently to mix for about 30 seconds.


• Add 3 drops of HI93703-51 Dispersing

• Insert the first cuvette (#1) with the reacted deionized water into the holder and ensure that the notch on the cap is positioned securely in the groove.

• Press Zero and the display will show the countdown prior to zeroing the blank. To skip the timer, press Zero twice. When the timer ends the meter will perform the zero reading. The display will show "-0.0-" when the meter is zeroed and ready for measurement.

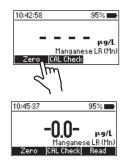




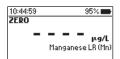


- Remove the cuvette.
- Insert the second cuvette (#2) with the reacted deionized water into the holder and ensure that the notch on the cap is positioned securely in the groove.



 Press Read to start the reading. The instrument displays the results in µg/L of manganese (Mn).

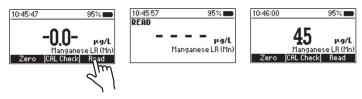

• Press the \blacksquare key and use the functional keys to select *Chemical Form*.

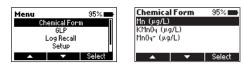

 Use the functional keys and press Select to change the displayed chemical formula to μg/L of potassium permanganate (KMnO₄) and permanganate (MnO₄⁻).



INTERFERENCES

Interference may be caused by:

- Calcium above 200 mg/L CaCO₃
- Magnesium above 100 mg/L $CaCO_3$
- Copper above 50 mg/L
- Nickel above 40 mg/L
- Aluminum, Cobalt above 20 mg/L
- Zinc above 15 mg/L
- Cadmium, Iron above 10 mg/L
- Lead above 0.5 mg/L

9. WARNING & ERROR DESCRIPTIONS

The instrument shows clear warning messages when erroneous conditions appear and when measured values are outside the expected range.

The information below provides an explanation of the errors and warnings, and recommended action to be taken.



Check the Zero cuvette

over limit. Wait for meter to cool down.

Continue

Meter temperature under limit. Put the meter in a warm place. Continue

Meter temperature

changing too fast. Redo Zero.

Continue

Warning Meter temperature

🗥 Warning

Warning


There is an excess amount of ambient light reaching the detector. Ensure that the notch on the cap is positioned securely in the groove before performing any measurements. If the issue persists, please contact Hanna Instruments technical support.

The sample and the zero cuvettes are inverted. Swap the cuvettes and repeat the measurement.

There is either too much light or the instrument can not adjust the light level. Please check the preparation of the zero cuvette and that the sample does not contain any debris.

The meter is either overheating or its temperature has dropped too low to operate within published accuracy specifications. The meter must be between 0 and 50 $^{\circ}$ C (32 and 122 $^{\circ}$ F) to perform any measurements.

Meter temperature has changed significantly since the zero measurement has been performed. The zero measurement must be performed again.

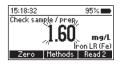

The measured value is outside the limits of the method. Verify that the sample does not contain any debris. Check the sample preparation and the measurement preparation.







Info	1
Tutorial Mode is Enabled.	
Continue	
	-

😧 Erroi	r
	rt the meter.
If iss	ue persists
contact te	echnical support.
	Continue

Date and time settings have been lost. Please reset the values. If the issue persists, please contact Hanna Instruments technical support.

English is the only available language. Help function is not available. Restart the meter. If the issue persists, please contact Hanna Instruments technical support.

Battery level is too low for the meter to function properly. Replace the batteries with new ones.


Tutorial mode has been enabled in the Setup menu. Press **Continue** and follow the prompt on the screen. Tutorial mode can be disabled in the Setup menu.

A critical error has occurred. Restart the meter. If the issue persists, please contact Hanna Instruments technical support.

10. BATTERY REPLACEMENT

To replace the instrument's batteries, follow these steps:

- Turn the instrument off by pressing and holding the 0 key.
- Remove the battery cover by turning it counterclockwise.
- Remove the old batteries, replace them with three new 1.5V AA batteries.
- Replace the battery cover, turn it clockwise to close.

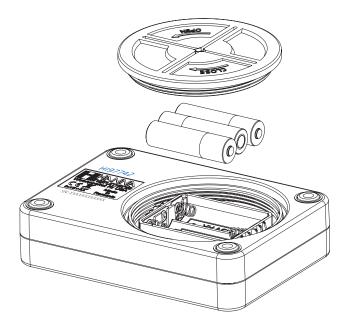

11. ACCESSORIES

11.1. REAGENT SETS

Code	Description
HI93746-01	Iron LR Reagent - 50 tests
HI93746-03	Iron LR Reagent - 150 tests
HI93748-01	Manganese LR Reagent - 50 tests
HI93748-03	Manganese LR Reagent - 150 tests

11.2. OTHER ACCESSORIES

Code	Description
HI7101413	blue carrying case for H1977xx and 3 CAL Check cuvettes
HI731318	cloth for wiping cuvettes (4 pcs.)
HI731331	glass cuvette (4 pcs.)
HI731336N	cap for glass cuvette (4 pcs.)
HI740143	1 mL graduated syringe (6 pcs.)
HI740144	pipette tip (6 pcs.)
HI740220	25 mL graduated glass vial (2 pcs.)
HI740229	100 mL graduated cylinder
HI93703-50	cuvette cleaning solution (230 mL)
HI97742-11	$\operatorname{CAL}\nolimits{\operatorname{Check}}^{\circledast}$ standards for Iron LR - cuvette kit
HI97748-11	$\operatorname{CAL}\nolimits{\operatorname{Check}}^{\circledast}$ standards for Manganese LR - cuvette kit

CERTIFICATION

All Hanna Instruments conform to the CE European Directives.



Disposal of Electrical & Electronic Equipment. The product should not be treated as household waste. Instead hand it over to the appropriate collection point for the recycling of electrical and electronic equipment which will conserve natural resources.

Disposal of waste batteries. This product contains batteries, do not dispose of them with other household waste. Hand them over to the appropriate collection point for recycling.

Ensuring proper product and battery disposal prevents potential negative consequences for the environment and human health. For more information, contact your city, your local household waste disposal service, the place of purchase or go to www.hannainst.com.



RECOMMENDATIONS FOR USERS

Before using this product, make sure it is entirely suitable for your specific application and for the environment in which it is used. Any variation introduced by the user to the supplied equipment may degrade the meter's performance. For yours and the meter's safety do not use or store the meter in hazardous environments.

WARRANTY

The H197742 is warranted for two years against defects in workmanship and materials when used for its intended purpose and maintained according to instructions. This warranty is limited to repair or replacement free of charge. Damage due to accidents, misuse, tampering or lack of prescribed maintenance is not covered.

If service is required, contact your local Hanna Instruments Office. If under warranty, report the model number, date of purchase, serial number (engraved on the bottom of the meter) and the nature of the problem. If the repair is not covered by the warranty, you will be notified of the charges incurred. If the meter is to be returned to Hanna Instruments, first obtain a Returned Goods Authorization (RGA) number from the Technical Service department and then send it with shipping costs prepaid. When shipping any meter, make sure it is properly packed for complete protection.

Hanna Instruments reserves the right to modify the design, construction or appearance of its products without advance notice.

World Headquarters

Hanna Instruments Inc. Highland Industrial Park 584 Park East Drive Woonsocket, RI 02895 USA www.hannainst.com



MAN97742

Printed in ROMANIA