Fig. 1

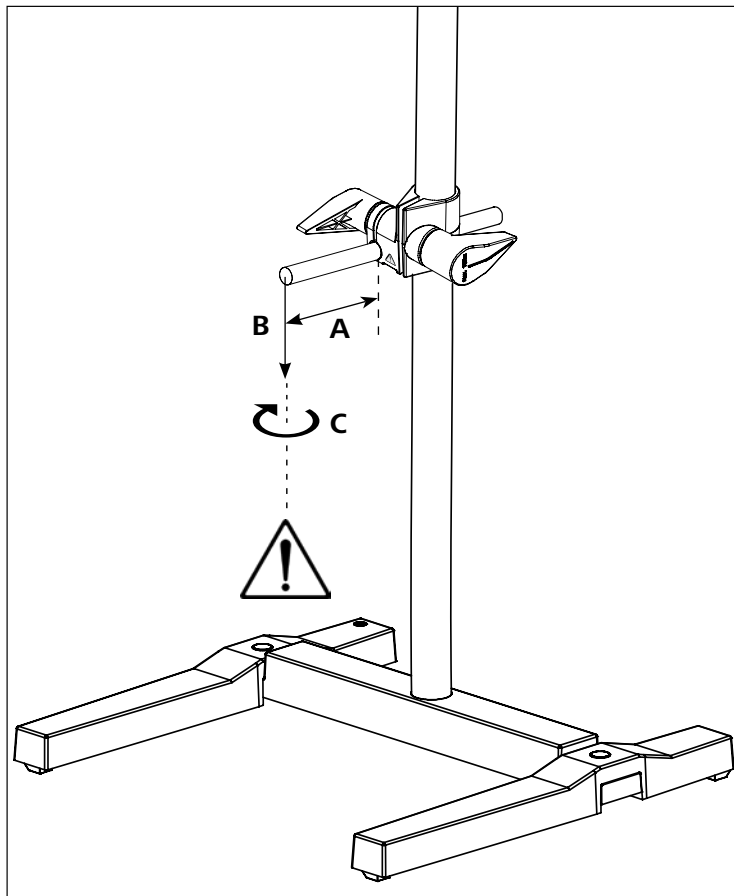
Fig. 2

IKA

designed for scientists

IKA-Werke GmbH & Co. KG
 Janke & Kunkel-Straße 10, 79219 Staufen, Germany
 Phone: +49 7633 831-0, Fax: +49 7633 831-98
 eMail: sales@ika.de

USA

IKA Works, Inc.
 Phone: +1 910 452-7059
 eMail: usa@ika.net

KOREA

IKA Korea Ltd.
 Phone: +82 2 2136 6800
 eMail: sales-lab@ika.kr

BRAZIL

IKA Brasil
 Phone: +55 19 3772 9600
 eMail: sales@ika.net.br

MALAYSIA

IKA Works (Asia) Sdn Bhd
 Phone: +60 3 6099-5666
 eMail: sales.lab@ika.my

CHINA

IKA Works Guangzhou
 Phone: +86 20 8222 6771
 eMail: info@ika.cn

POLAND

IKA Poland Sp. z o.o.
 Phone: +48 22 201 99 79
 eMail: sales.poland@ika.com

JAPAN

IKA Japan K.K.
 Phone: +81 6 6730 6781
 eMail: info_japan@ika.ne.jp

INDIA

IKA India Private Limited
 Phone: +91 80 26253 900
 eMail: info@ika.in

UNITED KINGDOM

IKA England LTD.
 Phone: +44 1865 986 162
 eMail: sales.england@ika.com

VIETNAM

IKA Vietnam Company Limited
 Phone: +84 28 38202142
 eMail: sales.lab-vietnam@ika.com

Discover and order the fascinating products of IKA online:
www.ika.com

IKAworldwide

IKAworldwide /// #lookattheblue

@IKAworldwide

Technical specifications may be changed without prior notice.

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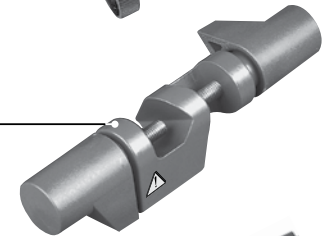
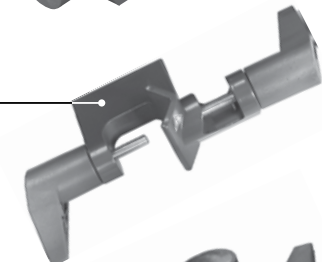
Safety instructions

- Read the mounting instructions in full before starting up and follow the safety instructions.
- Keep the operating instructions in a place where they can be accessed
- Safe operation of **IKA** laboratory devices with stands and cross sleeves is only guaranteed for combinations listed in the corresponding device operating instructions.
- Check that the laboratory device and the cross sleeve are properly tightened before each use.
- The laboratory device may only be moved on the stand when the device is at a standstill and unplugged!
- Clamped joints may loosen, resulting in danger to users and trial set-up!
- Hold the cross sleeve while loosening the clamped joints.
- While opening the clamped joints no sensitive parts are allowed to be under the device.
- Do not substitute turning handles or tighten with a tool.
- Do not use faulty cross sleeves!
- Observe the technical data!

Assembly

- Ensure that the extension arm on the laboratory device is tight; check regularly and retighten if necessary.
- Attach cross sleeve (**K**, see **Fig. 1**) to stand (**I**, see **Fig. 1**); open side of cross sleeve must face up (**R 182 / R 270**)!
- Mount laboratory device extension arm (**L**, see **Fig. 1**).
- When laboratory device is in desired position, firmly finger-tighten both turning handles (**M**, see **Fig. 1**) one after another.

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